



Instruction Manual

Growth and Maintenance of Mimic™ Insect Cells

Catalog no. 12552-014

Version B

21 August 2003

25-0540

Table of Contents

Kit Contents and Storage	iv
Accessory Products	1
Overview of Mimic™ Cells	2
Media.....	4
Cell Maintenance and Handling	5
Initiation of Cell Culture from Frozen Stock	8
Subculturing Cells in Adherent Culture	10
Subculturing Cells in Suspension Culture.....	11
Freezing Protocols	13
Support Protocols.....	14
Cell Log	16
Troubleshooting Guide.....	17
Purchaser Notification	22
Technical Service.....	23
References	25

Kit Contents and Storage

General Information

Mimic™ cells are prepared from Master Cell Bank cultures and shipped at less than 30 total passages.

Shipping and Storage

Shipping

Cells are shipped on dry ice.

Storage/Passaging

Store cells in liquid nitrogen. To initiate cell culture immediately upon receipt, see page 8.

Kit Contents

Catalog no. 12552-014 contains:

1 x 10⁷ cells in 1 ml 60% Grace's Media, 30% FBS, and 10% DMSO.

Product Qualification

Mimic™ insect cells are qualified using the following criteria:

- Cells must be in logarithmic growth with 98% viability before they are frozen
 - Cells must demonstrate that they can be recovered as healthy logarithmically growing cells within 2 to 3 days after thawing
 - Cells are tested independently and are certified to be free of mycoplasma
 - Expression of complex N-glycans containing galactose and sialic acid is verified by surface cell staining with FITC-conjugated lectins.
-

Accessory Products

Insect Culture Media

Mimic™ cells must be grown in Grace's Insect Medium, Supplemented, with serum. See below for ordering information:

Product	Amount	Catalog no.
Grace's Insect Medium, Supplemented	500 ml	11605-094
FBS, Heat-Inactivated	500 ml	10082-147
Grace's Insect Medium, Unsupplemented	500 ml	11595-030

Insect Expression Systems

Mimic™ cells can be used with a variety of insect expression systems; however, they are not compatible with Invitrogen's Bac-N-Blue™ Baculovirus Expression System (see note on page 3).

Product	Amount	Catalog no.
Bac-to-Bac® Expression System	1 kit	10359-016
InsectSelect™ Kit	1 kit	K800-01
InsectSelect™ BSD Kit	1 kit	K820-01

Transfection

For optimal results, we recommend using Cellfectin® Reagent for transfection.

Product	Amount	Catalog no.
Cellfectin® Reagent	1 ml	10362-010

Overview of Mimic™ Cells

Introduction

Mimic™ cells are transgenic insect cells that have been engineered to produce recombinant proteins with terminally sialylated N-glycans like those found in mammalian systems. These are identical to the SfSWT-1 cells developed by Donald L. Jarvis (Jarvis, *et al.*, 1998). This cell line is suitable for expressing recombinant glycoproteins with baculovirus and other insect expression systems, including the Bac-to-Bac® Expression System and the InsectSelect™ System.

This manual provides general information about the growth and maintenance of Mimic™ cell cultures.

Characteristics of Mimic Insect Cells

N-Glycosylation: Insect cells, like other eukaryotic cells, modify many of their proteins by N-glycosylation. However, endogenous insect cell N-glycan processing generally does not produce complex, terminally sialylated N-glycans like those produced by mammalian cells.

Mimic™ cell lines have been stably integrated with the following mammalian glycosyltransferases:

- α 2,6-sialyltransferase
- α 2,3-sialyltransferase (expressed, but not active)
- β 4-galactosyltransferase
- N-acetylglucosaminyltransferase I
- N-acetylglucosaminyltransferase II

These enzymes result in the production of highly processed, mammalian-like recombinant glycoproteins with biantennary, sialylated, complex glycans.

Origin: Mimic™ cells are derived from the Sf9 cell line, which are a traditional cell line used with baculovirus. Sf9 cells originated from the IPLBSF-21 cell line, derived from the pupal ovarian tissue of the fall army worm, *Spodoptera frugiperda* (O'Reilly *et al.*, 1992; Vaughn *et al.*, 1977).

Growth Characteristics: Mimic™ cells grow well in adherent and suspension culture. Media with serum is required. Doubling times for Mimic™ cells are similar to those for Sf9 cells.

Appearance: Mimic™ cells are spherical with some granular appearance.

Uses: Mimic™ cell lines are suitable for transfection and expression of recombinant proteins.

Insect Expression Systems

A number of insect expression systems are available from Invitrogen to facilitate expression of recombinant proteins in Mimic™ cells. The Bac-to-Bac® Baculovirus Expression System allows recombinant protein expression using baculovirus, while the InsectSelect™ System allows protein expression using a non-lytic, stable system. For more information about these expression systems, see our Web site (www.invitrogen.com) or call Technical Service (see page 22).

Continued on next page

Overview of Mimic™ Insect Cells, continued




Note

Mimic™ cells are not compatible with Invitrogen's Bac-N-Blue™ Baculovirus Expression System, because Bac-N-Blue™ viruses constitutively express active β -galactosidase, which will remove galactose from the N-glycan side chains of recombinant proteins expressed by Mimic™ cells.

N-Glycosylation Pathway

The figure below provides a model of insect cell and Mimic™ cell N-glycosylation pathways. Mimic™ cells have been modified to greatly enhance the production of complex, sialylated N-linked glycans, as shown in the lower right branch of the pathway.

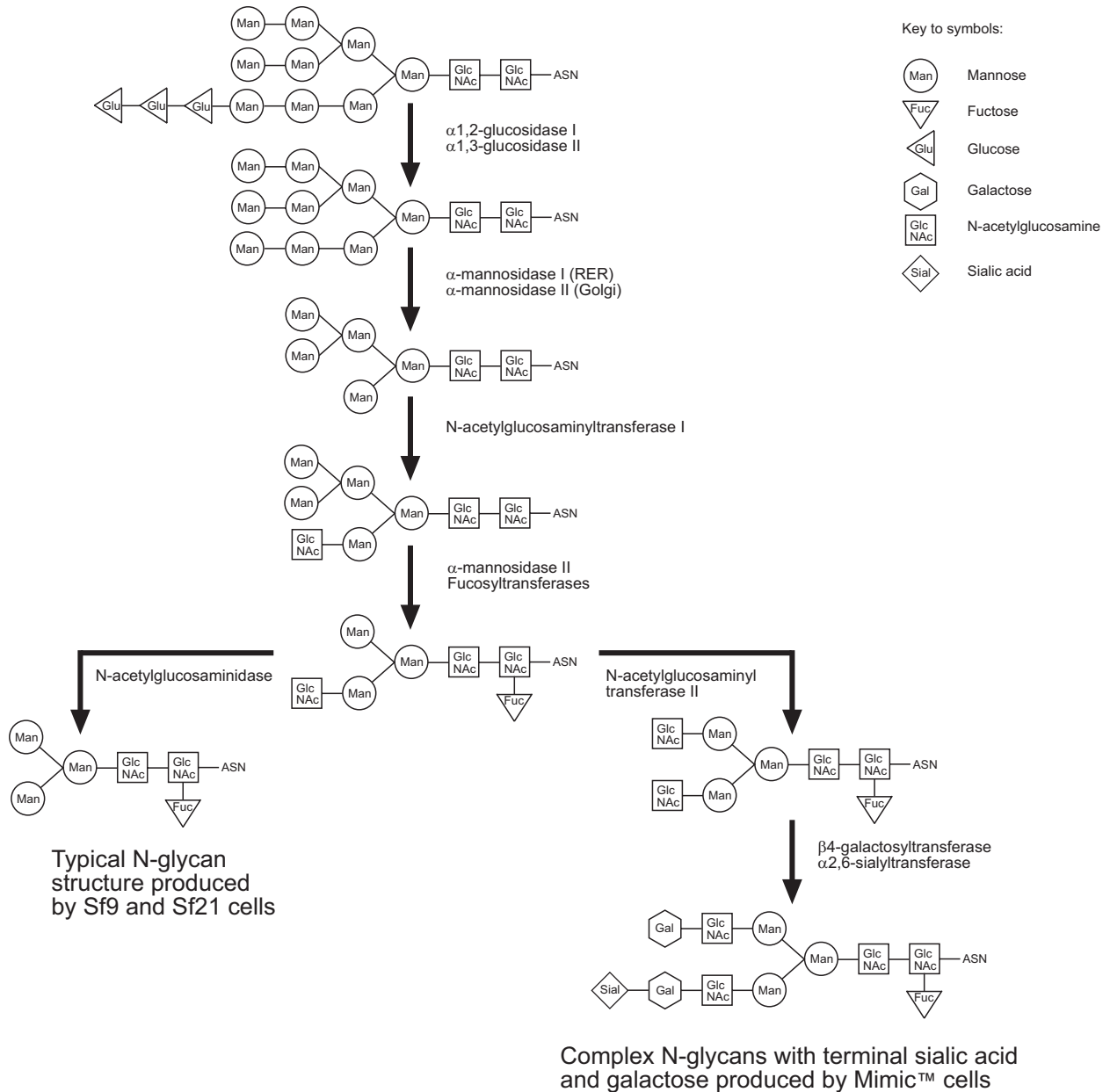

Media

Complete TNM-FH Mimic™ cells require complete TNM-FH (*Trichoplusia ni* Medium-Formulation Hink) for growth. TNM-FH is Grace's Insect Medium, Supplemented (available from Invitrogen—Cat. no. 11605-094), plus 10% fetal bovine serum (FBS, Cat. no. 10082-147). Serum provides a source of sialic acid and additional nutrients.

Grace's Insect Medium, Supplemented, includes TC yeastolate, lactalbumin hydrolysate, and L-glutamine.

When preparing complete TNM-FH, keep in mind the following:

- **Serum Sensitivity:** Serum does not need to be heat-inactivated prior to use. However, the quality of serum is important for optimal cell growth. We strongly recommend that you test a small aliquot of each new lot on your cells before using it. For more information on serum sensitivity, see **Troubleshooting Guide**, page 17.
 - **pH:** TNM-FH does not contain pH indicators. The normal pH for Mimic™ cells in this medium is 6.2. Unlike mammalian cell cultures, the pH rises gradually as the cells grow, but usually does not exceed pH 6.4.
 - **Stability:** Complete TNM-FH medium is stable for 1 month at +4°C.
-



Grace's Medium will not support the growth of Mimic™ cells without the additional supplements and serum.

Antibiotics

The use of antibiotics is optional. Many antibiotics are suitable for use with Mimic™ cells. The following table summarizes some of the most common.

Antibiotic	Working Conc.	Method of Action
Gentamycin	10 µg/ml	Inhibits bacterial protein synthesis
Amphotericin B (Fungizone)	0.25 µg/ml	Binds sterols and interferes with membrane permeability
Penicillin-Streptomycin	100-200 U/ml 100 µg/ml	Inhibits bacterial cell wall synthesis Inhibits bacterial protein synthesis

Preparation of Complete TNM-FH Medium

If possible, prepare complete TNM-FH medium just prior to use. The medium may be stored at +4°C for up to 1 month if it does not become contaminated.

1. Add 55 ml of FBS to 500 ml of Grace's Insect Medium, Supplemented. Add antibiotic, if desired, at the appropriate concentration.
 2. Filter this solution through a 0.2-µm filter into a sterile container.
 3. Store complete TNM-FH medium at +4°C.
 4. Equilibrate to room temperature before use.
-

Working With Baculovirus and Media

When working with recombinant or wild-type viral stocks, always maintain separate media bottles for cell culture and for virus work. Baculovirus particles can survive and be maintained in media at +4°C and will contaminate your cell cultures if added to plates or flask during passaging.

Cell Maintenance and Handling

Introduction	Use the recommended conditions, procedures, and techniques for subculturing Mimic™ cells.
Sterile Technique	All handling of Mimic™ cells should be performed in a laminar flow hood, using sterile conditions and equipment.
CO₂	CO ₂ exchange is not required for these cells.
Growth Temperature	<p>Fresh cell culture medium should be equilibrated to room temperature before use. Cells can be maintained at room temperature on the bench top or in a drawer; however, a 27°C controlled environment is recommended.</p> <p>Below 27°C, Mimic™ cells will grow more slowly, but will recover and resume normal doubling times once the temperature is returned to 27°C. Above 30°C, cells may display decreased viabilities. Cells under prolonged exposure to temperatures above 30°C should not be used.</p>
Confluency	Confluency is a marker for when to subculture your cells. A confluent monolayer is an adherent cell culture (dish, plate or flask) in which the cells have formed a single layer over the entire surface area available for growth.
Floaters	<p>Floaters are cells that are either loosely attached or suspended in the medium. Floaters are a normal occurrence, and are not necessarily an indicator of cell mortality.</p> <p>To check the viability of floaters, remove a small aliquot of medium containing floaters and assay for trypan blue exclusion (see protocol on page 15). If the viability is high (>95%), the cells may be propagated by transferring the medium containing floaters to a new, appropriately sized flask containing fresh medium.</p>
Determining Cell Density and Viability	<p>Follow the procedure below to determine total cell count and viability of cells in a culture:</p> <ol style="list-style-type: none">1. Transfer a small aliquot of the cell suspension culture to a microcentrifuge tube.2. Determine the cell density electronically using a Coulter Counter or manually using a hemacytometer chamber (see page 14).3. Determine cell viability using the trypan blue exclusion method (see page 15 for procedure). <p>Cell viability should be regularly assessed during cell counts in order to maintain optimal adherent and suspension cultures.</p>

Continued on next page

Cell Maintenance and Handling, continued

Minimum Viability Requirement

Cell viability should be at least 95% for healthy log-phase cultures. Cells below 95% viability are not growing under optimal conditions and should not be used in experiments. See **Troubleshooting Guide**, page 17 if your cell viability is below 95%.

Concentrating Cells

Cells can be concentrated if their density is too low (less than 5×10^5 cells/ml) to support log phase growth. Concentrating cells to a higher density (1×10^6 cells/ml) will induce log phase growth (see page 7).

Cell Log

We recommend keeping a record of cell density, viability, passage number, and other information as a decision-making and troubleshooting tool. See page 16 for a sample log form.

Removing and Adding Medium—Monolayers in Flasks

To **remove** medium from a flask of cells, tilt the flask so that all of the medium flows to one corner, away from the cell monolayer, and remove the medium carefully with a pipette. Avoid touching the cell monolayer.

To **add** medium to a flask: Gently pipet the medium down the side of the flask away from the cell monolayer.

Removing and Adding Medium—Monolayers in Plates

To **remove** medium from a plate containing a cell monolayer, tilt the plate at a 45°C angle so that the medium flows to one edge and aspirate the medium completely and very carefully using a Pasteur pipette.

To **add** medium to a plate, carefully and slowly add the medium against the side edge of the plate. Allow the medium to cover the entire surface of the plate, being careful not to dislodge cells.

Removing and Adding Medium—Suspension Cultures

Removing medium from a shake flask or spinner:

1. Remove the screw cap from one arm of the flask.
2. Carefully insert a pipette without touching the flask and remove medium from the culture.
3. Carefully remove pipette from the arm of the flask without touching the sides or dropping any medium.

Change pipettes each time one is inserted into the flask and then removed.

Adding medium to a shake flask or spinner:

1. Remove the screw cap from one arm of the flask.
2. Carefully insert pipette without touching the flask and add medium.
3. Carefully remove pipette from the arm of the flask without touching the sides or dropping any medium.

Change pipettes each time one is inserted into the flask and then removed.

Continued on next page

Cell Maintenance and Handling, continued

Concentrating Cells

To concentrate cells from a suspension culture (or resuspended cells from monolayer culture):

1. Transfer cell suspension to a sterile centrifuge tube of appropriate size and centrifuge for 5 minutes at $800 \times g$.
Important: Do not centrifuge at higher speeds. Mimic™ cells are very sensitive to centrifugal force.
 2. Carefully remove the supernatant without disturbing the pellet of cells.
 3. Add the desired volume of fresh, complete TNM-FH gently to the side of the tube and slowly pipette up and down 2 to 3 times to resuspend the cell pellet.
 4. Transfer to the desired, sterile container.
-

Cell Dispersion

Prior to performing transfections and plaque assays, cells need to be evenly distributed over the surface of a tissue culture plate. This ensures that:

- Cells do not distribute unevenly, leading to asymmetric monolayers.
- Maximum cell surface area is available for infection.

Procedure: To disperse cells, rock the flask or plate slowly by hand forward and backward, then side to side. Do this four times, watching carefully to be sure the liquid reaches all areas of the growth surface. Do not use a circular motion to disperse cells as this causes a concentration of cells around the edges of the plate rather than an even distribution

Initiation of Cell Culture from Frozen Stock

Introduction

Use the following protocols to initiate a culture from a frozen stock of Mimic™ cells.

Required Materials

- **Frozen Mimic™ cells**, in vials
 - **Water bath** at 37°C
 - **T-25 cm² flask** (per vial of cells)
 - **Complete TNM-FH** (see recipe on page 4)
 - **27°C constant-temperature incubator**
-

Sterile Technique

All handling of Mimic™ cells should be performed under sterile conditions in a laminar flow hood.

Experimental Outline

The table below outlines the steps to initiate a cell culture from frozen stock.

Step	Action
1	Remove cells from liquid nitrogen or dry ice (if just received)
2	Thaw cells into complete TNM-FH
3	Let cells attach for 30–45 minutes
4	Remove medium (with DMSO) and add fresh medium
5	Grow to confluency and passage
6	Subculture until cells are doubling every 24–30 hours and are 95% viable
7	Freeze down several vials of low passage cells as backup

Continued on next page

Initiation of Cell Culture from Frozen Stock, continued

Initiating Culture with Frozen Cells

Thawed Mimic™ cells do not always appear round; some may be amorphous or have a wrinkled appearance. A significant portion of the cells may have lysed. The cellular debris will be eliminated through successive rounds of subculturing.

The freezing process can damage cells, resulting in a 20–25% mortality rate upon thawing. For example, if 1×10^7 cells are frozen, $7.5\text{--}8 \times 10^6$ cells may be recovered.

1. Remove a vial of cells from dry ice or liquid nitrogen (if they were stored) and thaw rapidly in a 37°C water bath.
2. Remove from water bath and quickly decontaminate the outside of the vial by treating with 70% ethanol.
3. Transfer the 1 ml of cell suspension directly into a T-25 cm² flask containing 4 ml of complete TNM-FH.
4. Transfer flask to a 27°C incubator and allow cells to attach for 30–45 minutes.
5. After the cells are attached, gently remove the medium (see page 6). This must be done as soon as the cells have attached to remove the DMSO from the freezing medium. This step will also remove cellular debris and unhealthy cells that do not adhere.
6. Feed cells with 5 ml of fresh TNM-FH.
7. After 24 hours, change the medium. Continue to incubate until cells have formed a confluent monolayer.

At this point, cells should be subcultured in adherent culture (see following page). Mimic™ cells should be established in adherent culture before scaling up into suspension culture.

Subculturing Cells in Adherent Culture

Introduction

This section describes how to subculture Mimic™ cells in adherent culture. Mimic™ cells should be established in adherent culture before scaling up into suspension culture.

When to Subculture

Adherent cultures should be passaged at >80 percent confluency. **Do not overgrow cells.** Repeated subculturing of cells at densities past confluency will result in decreased doubling times and decreased viabilities.

We recommend thawing a fresh, low-passage culture of frozen cells every 3 months or 30 passages. We recommend keeping a log of the cell density, viability, passage number, and appearance. See page 16 for a sample cell log.

Required Materials

- T-flasks of desired size
 - Complete TNM-FH (see recipe on page 4)
 - Sterile pipettes
-

Recommended Dilution and Culture Volumes

Mimic™ cells should be diluted at no greater than a 1:2–1:3 dilution (volume of cells: final volume of medium) in order to maintain log phase growth.

Do not split cells back too far. Densities lower than 33% confluency will inhibit growth.

The table below summarizes recommended flask volumes for adherent culture.

Flask Size	Volume Range
T-25 cm ²	5–10 ml
T-75 cm ²	15–20 ml
T-150 cm ²	40–50 ml

Subculturing Procedure

We recommend sloughing to dislodge monolayers in adherent culture. Follow the procedure below to subculture adherent cultures:

1. Remove all but 5 ml of medium from flask (see page 6).
 2. Tilt flask on end so that the remaining medium flows to one corner, away from the cells.
 3. Draw up some of the remaining medium into a sterile Pasteur pipette and, starting at the bottom corner of the flask, stream medium across the cells (sloughing). Dislodge cells using side-to-side gentle streaming motion as you move from the bottom corner up to the top opposite corner of the flask.
 4. Determine cell density and viability (see protocols on pages 14–15).
 5. Seed cells at a 1:2–1:3 dilution in a T-flask (see sizes and volumes above) for adherent culture, or transfer to suspension culture at a density of 1×10^6 cells/ml (see following pages for information on suspension cultures).
-

Subculturing Cells in Suspension Culture

Introduction

This section describes how to establish and maintain Mimic™ cells in suspension culture.



Important

Mimic™ cells should be established in adherent culture before scaling up into suspension culture.

Required Equipment and Materials

For shaker cultures, you will need the following:

- **Adherent Mimic™ cells.** Cells must be grown to seed a shake flask at a minimum density of 1×10^6 cells per ml.
 - **27°C constant-temperature incubator;** CO₂ not required.
 - **Rotating, orbital shaker platform.**
 - **Disposable, sterile Erlenmeyer flasks.** Glass flasks without baffles may be used, but be sure to clean flasks thoroughly after each use to avoid potential toxicity.
 - **Complete TNM-FH containing Pluronic® F-68** at 0.1% final concentration.
-

Recommended Seeding Density and Culture Volumes

Suspension cultures should be seeded at a density of 1×10^6 cells/ml.

For shaker cultures, we recommend the flask sizes and total working volumes below:

<u>Flask Size</u>	<u>Volume Range</u>
125 ml	35–50 ml
250 ml	75–100 ml

Pluronic® F-68

For suspension cultures, Pluronic® F-68 (Catalog no. 24040-032) should be added to your complete TNM-FH medium at a final concentration of 0.1%.

Establishing Shake Flask Cultures

To initiate a suspension culture of Mimic™ cells in shake flasks:

1. Grow up enough log phase adherent cells to start a shake flask of desired size with 1×10^6 cells/ml. Determine cell density and viability of adherent cells according to the protocols on pages 14–15. Seeding at lower densities can cause decreased doubling times in the initial stages of culture.
Example: A 125-ml shake flask with 35 ml of culture at 1×10^6 cells/ml would require 3.5×10^7 viable cells.
 2. Seed a clean, sterile shake flask with cells at a density of 1×10^6 cells/ml. Use fresh, complete TNM-FH with 0.1% Pluronic®.
 3. Loosen flask caps to allow for aeration.
 4. Incubate flask at 27°C on a rotating orbital shaker at 85–95 rpm. Check cells daily for density and viability.
-

Subculturing Suspension Cell Cultures, continued

Subculturing Shake Flask Cultures

Cells in suspension culture should be passaged when they reach a density of 2.0 to 2.5×10^6 cells/ml.

Example: A 125-ml shake flask containing 50 ml of cell culture is at a density of 2×10^6 cells/ml. Remove 25 ml of medium and cells and seed in a new flask with 25 ml of fresh complete medium for a final density of 1×10^6 cells/ml.

Maintaining Shake Flask Cultures

- **Renew cultures every month:** Cells can be kept in suspension for up to one month in log phase. At this point, a new culture should be initiated as cells begin to lose their infectivity and are no longer optimal.
 - **Maintain log phase cell density:** Maintain density above 1×10^6 cells/ml for log phase growth. When cells reach a density of 2.0 to 2.5×10^6 cells/ml, they should be diluted with fresh medium to 1×10^6 cells/ml.
 - **Maintain volumes for adequate aeration:** The total culture volume in a shake flask should not exceed 40% of the indicated volume of the flask for proper aeration (e.g., a 250-ml flask should not contain more than 100 ml of culture).
 - **To reduce accumulation of cell debris and metabolic waste by-products,** we recommend gently centrifuging the cell suspension at $100 \times g$ for 5 minutes and resuspending the cell pellet in fresh, complete TNM-FH once every 3 weeks (see page 7 for detailed instructions on concentrating cells).
 - **Keep a log:** We recommend keeping a log of the cell density, viability, passage number, and appearance. See page 16 for a sample cell log.
-

Spinner Cultures

You can scale up from shake flasks to spinner cultures using the guidelines below.

- **Flask:** We recommend starting with 100-ml or 250-ml spinner flasks, as these require fewer total cells than larger flasks to initiate culture.
 - **Seeding density:** 1×10^6 cells per ml.
 - **Spinner culture volume:** To ensure proper aeration:
 - The **maximum** culture volume should not exceed 50% of the spinner volume (e.g., a 250-ml spinner should have a culture volume of ≤ 125 ml).
 - The **minimum** culture volume should submerge the impeller to a depth of at least 1 cm (e.g., a 250-ml spinner should have a culture volume of ≥ 80 ml).
 - **Impeller:** We recommend using a spinner flask with a **vertical impeller** rather than one with a hanging stir-bar assembly, as the vertical impeller provides better aeration. Incubate at 27°C with an impeller speed of 80–95 rpm.
 - **Spinner Caps:** Finger-tighten caps. Loose caps are not necessary for aeration.
 - **Subculturing:** Subculture cells when density reaches 2.0 to 2.5×10^6 cells/ml. Maintain density above 1×10^6 cells/ml for log phase growth.
 - **Cleaning:** Do not use soap when cleaning spinners, as residue will cause cell mortality. See page 15 for more information.
-

Freezing Protocols

Introduction

Once Mimic™ cell lines are established and doubling regularly, they can be frozen. Cells should be at least 90% viable and in mid-log exponential growth. You will need enough cells to freeze down 2–4 cryovials at a density of 1×10^7 each. We recommend that you freeze down several vials as soon as your cell culture meets these requirements (e.g., at as low a passage number as possible).

Freezing Procedure

To make the freezing medium, combine:

60% Grace's Insect Medium, Unsupplemented

30% FBS

10% DMSO

Note that Grace's Insect Medium does not contain supplements or FBS.

1. Count cells using a hemacytometer, as described on page 14. You need enough cells to freeze down 2–4 cryovials at a density of 1×10^7 each. Cells can come from suspension or adherent culture.
2. Set up sterile cryovials on ice. Be sure to label the cryovials.
3. Centrifuge the cells at $400\text{--}600 \times g$ for 10 minutes at room temperature. Remove the supernatant.
4. Resuspend cells to a density of 1×10^7 in the freezing medium.
5. Transfer 1 ml of the cell suspension to sterile cryovials.
6. Freeze cells in an automated, controlled-rate freezing apparatus or using a manual method following standard procedures. For ideal cryopreservation, the temperature should decrease at a rate of 1°C per minute.
7. Store in liquid nitrogen.

Check the viability and recovery of frozen cells 24 hours after storing vials in liquid nitrogen by following the procedure on page 8 for initiating culture from frozen stock.

Support Protocols


Introduction

Guidelines to maintain and optimize Mimic™ cell cultures are provided in this section. Included are protocols for cleaning spinner flasks.

Determining Cell Density— Hemocytometer

Hemocytometers may be obtained from most major laboratory suppliers (e.g., Baxter Scientific). The procedure below provides some general directions on how to use the hemocytometer. For more details, refer to King and Possee, 1992.

1. Clean the chamber and coverslip with alcohol. Dry and fix the coverslip in position.
2. Harvest cells. Add 10 μl of the cells to the hemocytometer. Do not overflow.
3. Place chamber in the inverted microscope under the 10X objective. Use phase contrast to distinguish cells.
4. Count the cells in the large, central gridded square (1 mm^2). The gridded square is circled in the graphic below. Multiply by 10^4 to estimate the number of cells per ml. Prepare duplicate samples and average the count.

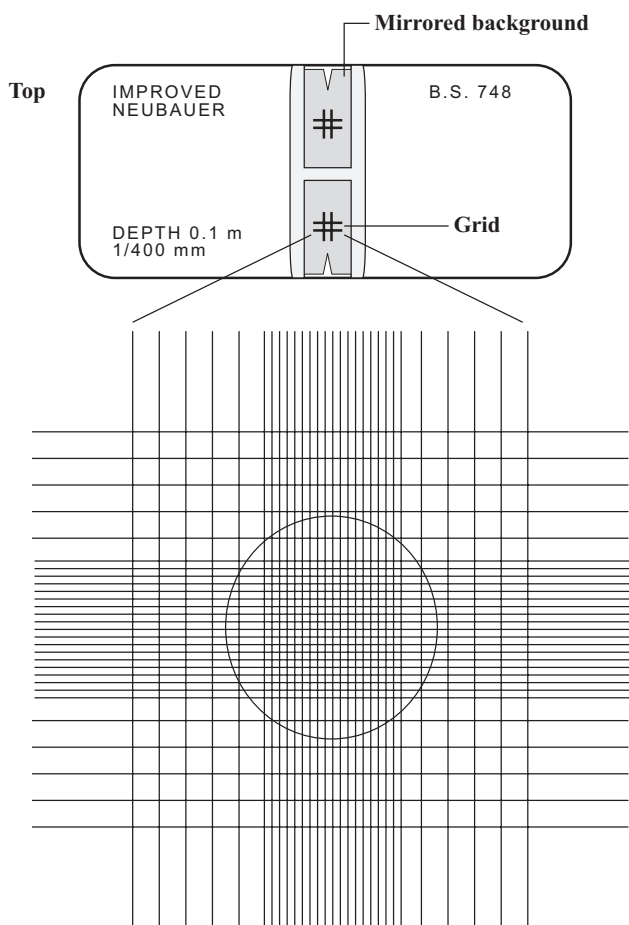

Continued on next page

Support Protocols, continued

Determining Cell Viability—Trypan Blue Exclusion

Trypan blue dye molecules are excluded by the membranes of viable cells, but readily enter dead cells. After staining with trypan blue, cells that are blue are considered nonviable.

The following procedure will enable you to quickly and accurately determine cell viability. Calculate viability by dividing the number of viable (non-blue) cells by the total number of cells in the grids on a hemacytometer.

1. Determine the cell density of your cell suspension using a hemacytometer (see previous page).
2. Prepare a 0.4% solution of trypan blue in buffered isotonic salt solution, pH 7.2 to 7.3 (i.e., phosphate-buffered saline).
3. Add 0.1 ml of trypan blue stock to 1 ml of cells.
4. Load a hemacytometer and examine immediately under a microscope at low magnification.
5. Count the number of blue staining cells and the number of total cells. Cell viability should be at least 95% for healthy log-phase cultures.

$$\% \text{ viable cells} = \left[1.00 - \left(\frac{\text{Number of blue cells}}{\text{Number of total cells}} \right) \right] \times 100$$

To calculate the number of viable cells per ml of culture, use the formula below. Remember to correct for the dilution factor.

$$\text{Number of viable cells} \times 10^4 \times 1.1 = \text{cells/ml culture}$$

Cleaning Spinner Flasks

Proper spinner care is important for both cell health and to prevent cross-contamination of your cultures with different viral stocks. Baculovirus particles can survive a cycle through the autoclave. We have developed a protocol that uses two cycles in the autoclave to eliminate virus contamination of future spinner cultures.

1. Wash with 7X glass cleaner (Bellco), or substitute 10% acetic acid if necessary. Wash while spinning, approximately 2 hours. The 7X glass cleaner is preferable, as trace amounts are not detrimental to the cells. Acetic acid can be detrimental if not rinsed away completely. **Do not use detergents.**
 2. Rinse 5 times with tap water.
 3. Rinse 5 times with deionized water.
 4. Tighten spinner cap and o-ring around impeller bar; spinner arm caps should be finger-tightened.
 5. Autoclave once wet with deionized water for 45 minutes on liquid cycle.
Note: You can repeat this step more than once if desired.
 6. Autoclave once dry for 45 minutes on dry cycle.
-

Cell Log

Cell Line: _____ Cell Lot #/Date of Freeze down: _____

Date of Initiation: _____ Medium: _____ Medium Lot No.: _____

Passage No.	Date of Passage	Cell Density	Cell Viability	Comments
1				
2				
3				
4				
5				
6				
7				
8				
9				
10				
11				
12				
13				
14				
15				
16				
17				
18				
19				
20				
21				
22				
23				
24				
25				
26				
27				
28				
29				
30				

Record any changes in medium and medium lot number in the "Comment" column as well as any notes about cell appearance.

Troubleshooting Guide

Introduction

The following section summarizes the most common problems associated with insect cell culture. Morphological changes in the cells or changes in the growth rate can indicate an underlying problem with the culture. If you have kept a record of cell viabilities and doubling times, this can be helpful in diagnosing a problem when it comes up. Use this table as a general guide for potential problems—what might cause them and what to do to solve and prevent them.

Adherent Cells: Morphology Changes—First Week in Culture

Problem	Potential Causes	Solutions
Cells are granular and/or floating	Mimic™ cells may become loosely attached or float, even when viable.	Transfer cells (including floaters) to suspension culture after required density has been attained (1×10^6 cells/ml). Check viability of floaters using trypan blue exclusion method (page 15).
	Medium was not removed within 1 hour of thawing cells. DMSO in the freezing medium can be harmful to cells.	Remove medium, add fresh medium.
	Cells are exhibiting serum sensitivity.	Try a new type or lot of serum, thaw new cells.
Cell lysis or debris	Aging cell culture with passage number greater than 30.	Thaw new cells of lower passage number.

Continued on next page

Troubleshooting Guide, continued

Adherent Cells: Morphology Changes—More Than One Week in Culture

Problem	Potential Causes	Solutions
Decreased viability and growth rate	Contamination with wild-type or recombinant baculovirus. You will see lysis, debris and/or swelling of cells. Inclusions in the nucleus indicates wild-type infection.	Discard the old culture and initiate a fresh culture. Thaw new cells. Decontaminate hood and equipment. Make up new medium.
	Too much mechanical manipulation during subculturing.	Thaw new cells. Switch to another method of subculturing (e.g., sloughing).
Floater (greater than 5-10% of cells in culture)	Mimic™ cells may become loosely attached or float, even when viable.	Transfer cells (including floaters) to suspension culture after required density has been attained (1×10^6 cells/ml). Check viability of floaters using trypan blue exclusion method (page 15).
	Passaging too often before confluency.	Remove medium / floaters. Replace with fresh medium. Passage only at confluency.
	Cells are overgrown.	If this is the first time this has occurred, replace medium and split cells. Thaw new cells if this has happened more than once.
	Sensitivity to a new lot or brand of serum (FBS).	Test your culture with another serum. We recommend using serum from GIBCO®.
Cells swollen, spots in the nucleus	Contamination with wild-type or recombinant baculovirus.	Discard the old culture and initiate a fresh culture. Thaw new cells. Make new medium. Decontaminate equipment.

Continued on next page

Troubleshooting Guide, continued

Adherent Cells: Growth and/or Viability Decrease

Problem	Potential Causes	Solutions
Cell doubling time is too long (>30 hours)	Cells have grown past confluency more than once.	Thaw new cells and check your cell cultures daily to avoid overgrowth.
	Cells are at a high passage number (greater than 30 passages).	Thaw new cells and make sure that you are freezing down cells at a low passage number (less than 10) so that new cultures are optimal.
	Cells are repeatedly being passaged before confluency.	Let cells grow to confluency for next passage. If doubling time does not increase, thaw new cells.
	Cells are split back below 20% confluency repeatedly.	Concentrate cells to a confluency of 50% or greater and plate out (see page 7).
Cell viability is less than 90%	Subculturing technique too harsh.	Thaw new cells and try using a different subculturing technique (e.g., sloughing, page 10).
	Bacterial or fungal contamination.	Discard the old culture and initiate a fresh culture. Decontaminate equipment. Prepare fresh medium.
	Contamination with wild-type or recombinant baculovirus. You will see lysis, debris and/or swelling of cells. Inclusions in the nucleus indicates wild-type infection.	Discard the old culture and initiate a fresh culture. Decontaminate equipment. Prepare fresh medium. Keep separate bottles of medium for cell culture and for virus work. Do not work with cell culture at the same time that you are working with virus in the tissue culture hood.

Continued on next page

Troubleshooting Guide, continued

Suspension Culture: Morphology Changes

Problem	Potential Causes	Solutions
Clumping of cells	Shake/spin rate is too slow.	Use 80–95 rpm.
Cell lysis or debris	Aging cell culture has been maintained in suspension longer than two months.	Thaw new cells of low passage number and initiate a new suspension culture.
	Cell shearing due to shaking/spinning.	Add Pluronic® F-68 at 0.1% final concentration.
	Contamination with wild-type or recombinant baculovirus. You will see lysis, debris and/or swelling of cells. Inclusions in the nucleus indicates wild-type infection.	Discard the old culture and initiate a fresh culture. Thaw new cells and start a new culture. Decontaminate hood and equipment. Make up new medium. Keep separate bottles of medium for cell culture and viral work.
Cells swollen, spots in the nucleus	Contamination with wild-type or recombinant baculovirus.	See above recommendations for baculovirus contamination.

Suspension Culture: Growth and/or Viability Decrease

Problem	Potential Causes	Solutions
Cell doubling time is too long (>30 hours)	Culture seeded or split back to a cell density that is too low (less than 5×10^5 cell/ml).	Concentrate cells to 1×10^6 cell/ml (see page 7). This will boost them into log phase growth.
	Spinner overgrown to a cell density $>3 \times 10^6$ cells/ml.	<ol style="list-style-type: none"> 1. Split cells to a density of 1.5×10^6 cells/ml. 2. Grow them overnight to a density of 2.0 to 2.5×10^6 cells/ml. 3. Split them back to a density of 1.0×10^6 cells/ml, continue normal maintenance (see page 15).

Continued on next page

Troubleshooting Guide, continued

Spinner Culture: Growth and/or Viability Decrease

Problem	Potential Causes	Solutions
Cell doubling time is too long (>30 hours), continued	Inadequate aeration.	The volume of culture should not exceed 40% the designated capacity of the shake flask, or 50% the capacity of the spinner.
		Add Pluronic® F-68 at 0.1% final concentration.
		Volume of culture should meet minimum volume requirements (i.e., 1×10^6 cells/ml).
		For spinner cultures, impeller should spin with a smooth continuous motion. Jerky or jumping motion of the impeller does not provide good aeration.
Viability < 90%	Cell shearing due to shaking/spinning.	Shaker/spinner speed is too high; recommended speed is 80–95 rpm. Add Pluronic® F-68 at 0.1% final concentration.
	Inadequate aeration.	See solutions above for Inadequate Aeration .
	Bacterial or fungal Contamination.	Discard the old culture and initiate a fresh culture. Decontaminate hood and equipment. Make up new medium. Keep separate bottles of medium for cell culture and viral work. Try adding antibiotics to your medium.
	Contamination wild-type or recombinant baculovirus. You will see lysis, debris and/or swelling of cells. Inclusions in the nucleus indicates wild-type infection.	Discard the old culture and initiate a fresh culture. Adequate sterilization of spinner flasks between uses where there is contact with baculovirus is required. See page 15.

Purchaser Notification

Information for European Customers

The Mimic™ cell line is genetically modified and contains the following stably integrated, pUC-derived plasmids: pIE1Hygro, pIE1GlcNAcTI, pIE1GlcNAcTII, pIE1ST6GalI, and pIE1ST3GalIV (Hollister, *et al*). As a condition of sale, this product must be in accordance with all applicable local legislation and guidelines including EC Directive 90/219/EEC on the contained use of genetically modified organisms.

Limited Use Label License No. 58: Stably Transformed Insect Cell Lines and Baculovirus Vectors

This product is the subject of one or more of U.S. Patent Nos. 5,077,214 and 5,162,222 licensed to Invitrogen Corporation by The Texas A & M University System. This product is sold for research purposes only. A separate license under the rights of The Texas A & M University system is required for use of this product for other than for research purposes. Inquiries should be directed to: Technology Licensing Office, Texas A & M University, MS 3369, College Station, TX 77843. Tel: 409-847-8682; Fax: 409-845-1402.

Limited Use Label License No. 139: Glycosylating Cell Lines

This product is the subject of U.S. Patent No. 6,461,863 assigned to University of Wyoming and licensed exclusively to Invitrogen. The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) not to transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. Invitrogen Corporation will not assert a claim against the buyer of infringement of the above patent(s) based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is willing to accept return of the product with a full refund. For information on purchasing a license to this product for purposes other than research, contact Licensing Department, Invitrogen Corporation, 1600 Faraday Avenue, Carlsbad, California 92008. Phone (760) 603-7200. Fax (760) 602-6500.

Technical Service

World Wide Web



Visit the Invitrogen Web Resource using your World Wide Web browser. At the site, you can:

- Get the scoop on our hot new products and special product offers
- View and download vector maps and sequences
- Download manuals in Adobe® Acrobat® (PDF) format
- Explore our catalog with full color graphics
- Obtain citations for Invitrogen products
- Request catalog and product literature

Once connected to the Internet, launch your Web browser (Internet Explorer 5.0 or newer or Netscape 4.0 or newer), then enter the following location (or URL):

<http://www.invitrogen.com>

...and the program will connect directly. Click on underlined text or outlined graphics to explore. Don't forget to put a bookmark at our site for easy reference!

Contact Us

For more information or technical assistance, call, write, fax, or email. Additional international offices are listed on our Web page (www.invitrogen.com).

Corporate Headquarters:

Invitrogen Corporation
1600 Faraday Avenue
Carlsbad, CA 92008 USA
Tel: 1 760 603 7200
Tel (Toll Free): 1 800 955 6288
Fax: 1 760 602 6500
E-mail:
tech_service@invitrogen.com

Japanese Headquarters:

Invitrogen Japan K.K.
Nihonbashi Hama-Cho Park Bldg. 4F
2-35-4, Hama-Cho, Nihonbashi
Tel: 81 3 3663 7972
Fax: 81 3 3663 8242
E-mail: jpinfo@invitrogen.com

European Headquarters:

Invitrogen Ltd
Inchinnan Business Park
3 Fountain Drive
Paisley PA4 9RF, UK
Tel: +44 (0) 141 814 6100
Tech Fax: +44 (0) 141 814 6117
E-mail: eurotech@invitrogen.com

MSDS Requests

To request an MSDS, visit our Web site at www.invitrogen.com. On the home page, go to 'Technical Resources', select 'MSDS', and follow instructions on the page.

Continued on next page

Technical Service, continued

Limited Warranty

Invitrogen is committed to providing our customers with high-quality goods and services. Our goal is to ensure that every customer is 100% satisfied with our products and our service. If you should have any questions or concerns about an Invitrogen product or service, contact our Technical Service Representatives.

Invitrogen warrants that all of its products will perform according to specifications. The company will replace, free of charge, any product that does not meet those specifications. This warranty limits Invitrogen Corporation's liability only to the cost of the product. No warranty is granted for products beyond their listed expiration date. No warranty is applicable unless all product components are stored in accordance with instructions. Invitrogen reserves the right to select the method(s) used to analyze a product unless Invitrogen agrees to a specified method in writing prior to acceptance of the order.

Invitrogen makes every effort to ensure the accuracy of its publications, but realizes that the occasional typographical or other error is inevitable. Therefore Invitrogen makes no warranty of any kind regarding the contents of any publications or documentation. If you discover an error in any of our publications, please report it to our Technical Service Representatives.

Invitrogen assumes no responsibility or liability for any special, incidental, indirect or consequential loss or damage whatsoever. The above limited warranty is sole and exclusive. No other warranty is made, whether expressed or implied, including any warranty of merchantability or fitness for a particular purpose.

References

- Davis, T. R., Trotter, K. M., Granados, R. R., and Wood, H. A. (1992) Baculovirus Expression of Alkaline Phosphatase as a Reporter Gene for Evaluation of Production, Glycosylation, and Secretion. *Bio/Technology* 10, 1148-1150.
- Granados, R. R., Guoxun, L., Derksen, A. C. G., and McKenna, K. A. (1994) A New Insect Cell Line from *Trichoplusia ni* (BTI-Tn-5B1-4) Susceptible to *Trichoplusia ni* Single Enveloped Nuclear Polyhedrosis Virus. *J. Invertebr. Pathol.* 64, 260-266.
- Hink, W. F., Thomsen, D. R., Davidson, D. J., Meyer, A. L., and Castellano, F. J. (1991) Expression of Three Recombinant Proteins using Baculovirus Vectors in 23 Insect Cell Lines. *Biotechnol. Prog.* 7, 9-14.
- Hollister, J.R., Grabenhorst, E., Nimtz, M., Conradt, H., and Jarvis, D.L. Engineering the protein N-glycosylation pathway in insect cells for production of biantennary, complex N-glycans. *Biochemistry*. *In press*.
- Jarvis, D.L. and Finn, E.F. (1996) Modifying the insect cell N-glycosylation pathway with immediate early baculovirus expression vectors. *Nat. Biotechnol.* 14, 1288-1292.
- Jarvis, D.L., Fleming, J.G.W., Kovacs, G.R., Summers, M.D. and Guarino, L.A. (1990) Use of early baculovirus promoters for continuous expression and efficient processing of foreign gene products in stably transformed lepidopteran cells. *Bio/Technology* 8, 950-955.
- Jarvis, D.L., Kowar, Z.S., and Hollister, J.R. (1998) Engineering N-glycosylation pathways in the baculovirus-insect cell system. *Curr. Opin. Biotechnol.* 9: 528-33
- King, L. A., and Possee, R. D. (1992) *The Baculovirus Expression System: A Laboratory Guide* (New York, NY: Chapman and Hall).
- O'Reilly, D. R., Miller, L. K., and Luckow, V. A. (1992) *Baculovirus Expression Vectors: A Laboratory Manual* (New York, N. Y.: W. H. Freeman and Company).
- Richardson, C. D. (1995) Baculovirus Expression Protocols. In *Methods in Molecular Biology*, Vol. 39. (J. M. Walker, ed.) Humana Press, Totowa, NJ.
- Seo, N.S., Hollister, J.R., and Jarvis, D.L. (2001) Mammalian glycosyltransferase expression allows sialoglycoprotein production by baculovirus-infected insect cells. *Protein Expr. Purif.* 22: 234-41
- Vaughn, J. L., Goodwin, R. H., Tompkins, G. J., and McCawley, P. (1977) The Establishment of Two Cell Lines from the Insect *Spodoptera frugiperda* (Lepidoptera: Noctuidae). *In Vitro* 13, 213-217.
- Wickham, T. J., Davis, T., Granados, R. R., Shuler, M. L., and Wood, H. A. (1992) Screening of Insect Cell Lines for the Production of Recombinant Proteins and Infectious Virus in the Baculovirus Expression System. *Biotechnol. Prog.* 8, 391-396.
- Wickham, T. J., and Nemerow, G. R. (1993) Optimization of Growth Methods and Recombinant Protein Production in BTI Tn-5B1-4 Insect Cells using the Baculovirus Expression Vector. *Biotechnol. Prog.* 9, 25-30.

©2002-2003 Invitrogen Corporation. All rights reserved.

For research use only. Not intended for any animal or human therapeutic or diagnostic use.



United States Headquarters:

Invitrogen Corporation
1600 Faraday Avenue
Carlsbad, California 92008
Tel: 1 760 603 7200
Tel (Toll Free): 1 800 955 6288
Fax: 1 760 603 7229
Email: tech_service@invitrogen.com

European Headquarters:

Invitrogen Ltd
3 Fountain Drive
Inchinnan Business Park
Paisley PA4 9RF, UK
Tel (Free Phone Orders): 0800 269 210
Tel (General Enquiries): 0800 5345 5345
Fax: +44 (0) 141 814 6287
Email: eurotech@invitrogen.com

International Offices:

Argentina 5411 4556 0844
Australia 1 800 331 627
Austria 0800 20 1087
Belgium 0800 14894
Brazil 0800 11 0575
Canada 800 263 6236
China 10 6849 2578
Denmark 80 30 17 40

France 0800 23 20 79
Germany 0800 083 0902
Hong Kong 2407 8450
India 11 577 3282
Italy 02 98 22 201
Japan 03 3663 7974
The Netherlands 0800 099 3310
New Zealand 0800 600 200
Norway 00800 5456 5456

Spain & Portugal 900 181 461
Sweden 020 26 34 52
Switzerland 0800 848 800
Taiwan 2 2651 6156
UK 0800 838 380

For other countries see our website

www.invitrogen.com

