

Protein Sample Preparation and Analysis Application Manual



Pall is the world leader in the design, production and supply of membranes, filtration devices, and separation systems. Pall's cutting edge filtration technologies make it possible to manufacture more effective medicines, better-tasting food, more efficient electronic devices, and much more. Wherever there is a need for a pure product, there is a need for Pall's technologies.

From multiple manufacturing facilities located in North America, Europe, and Asia, Pall provides enabling and process enhancing technologies to industries ranging from medical/pharmaceutical, aerospace, automotive, environmental, microelectronics, and semiconductors, to municipal and industrial water treatment, fuels, chemicals, energy, and food and beverages. Pall holds an impressive portfolio of products, quality systems, and experience to meet the diverse and demanding requirements of our customers. Our products are a direct result of consideration of your needs and smart collaborations.

Pall Life Sciences partners with our customers in the search for solutions to real-world health issues. Our vast product portfolio, practical knowledge, and technical know-how enable us to deliver products and services that truly make a difference. From innovations in drug discovery processes to ensuring safe water supplies, Pall technologies touch people's lives.

Contact Information

Pall Life Sciences separation and detection products can be ordered through Pall offices and distributors worldwide. For information contact:

PALL Life Sciences

**600 South Wagner Road
Ann Arbor, MI 48103-9019 USA**

**Tel: 800.521.1520 (in USA)
734.665.0651 (global)**

Fax: 734.913.6114

Customer Service:

**Tel: 800.521.1520 (in USA)
734.913.6194 (global)**

Fax: 734.913.6495

Technical Service:

**Tel: 800.521.1520 (in USA)
734.913.6123 (global)**

Fax: 734.913.6576

Email: Lab@pall.com

www.pall.com/lab

1.0 INTRODUCTION	
1.1 Proteomics Process Flow	1
1.2 Product Selection	2
2.0 PRE-ANALYTICAL APPLICATIONS	
2.1 Abundant Protein Removal	7
2.1.1 Introduction	7
2.1.2 Enchant™ Albumin Depletion Kit	10
2.1.3 Enchant IgG Purification Kit	19
2.1.4 Enchant Multi-Protein Affinity Separation Kit	33
2.1.5 BioSeptra® Chromatography Resins	43
2.2 Protein Fractionation (Complexity Reduction)	47
2.2.1 Introduction	47
2.2.2 Q, S, DEAE, and CM Ion Exchange Media	49
2.2.2.1 Ceramic HyperD® Ion Exchange Resin	49
2.2.2.2 Mustang® Ion Exchange Membrane	65
2.2.3 Affinity Separations – Chromatography Resins	72
2.2.3.1 Heparin HyperD M Resin	72
2.2.3.2 Lysine HyperD Resin	80
2.2.3.3 IMAC HyperCel™ Resin for Immobilized Metal Ion Affinity Chromatography	88
2.2.3.4 Protein A Ceramic HyperD F Resin	100
2.2.4 Mixed Mode	112
2.2.4.1 MEP HyperCel Mixed Mode Resin	112
2.2.4.2 HA Ultrogel® Resin	121
2.2.5 Size Fractionation	132
2.2.5.1 Ultrogel AcA Size Exclusion Chromatography Resin	132
2.3 Detergent Removal	141
2.3.1 SDR HyperD Solvent and Detergent Removal Chromatography Resin	141
2.4 Concentration, Desalting, and Buffer Exchange	152
2.4.1 Introduction	152
2.4.2 Concentration and Desalting Using Ultrafiltration	154
2.4.2.1 Concentration and Diafiltration of Samples (< 60 mL) in a Spin Filter	154
2.4.2.2 Desalting and Buffer Exchange of Samples (< 1 mL) in an AcroPrep™ 96 and 384 Multi-Well Filter Plate	179
2.4.3 Desalting with Size Exclusion Media	188
2.4.3.1 Ultrogel AcA 202 and Trisacryl® GF-05M Resins	188
2.5 Generic Clarification of Samples by Microporous Filtration (Particulate Removal)	198
2.5.1 Introduction	198
2.5.2 AcroPrep 96 and 384 Multi-Well Filter Plate for 0.1-1.0 mL Samples	200
2.5.3 Clarification of Samples (< 1 mL) in Spin Filters (Nanosep® Centrifugal Devices)	210
2.5.4 Clarification of Samples (1-100 mL) in an Acrodisc® Syringe Filter	214
2.5.5 Filtration of Reagents (> 50 mL) with a VacuCap® Vacuum Filtration Device	221
2.5.6 Clarification of Plasma or Serum Samples to Remove Cryoprecipitate Using Acrodisc Syringe Filters with a Prefilter	226
2.5.7 Prefiltration of Plasma or Serum Samples Using Acrodisc Syringe Filters with Glass Fiber	229
2.6 Endotoxin Removal	231
2.6.1 Endotoxin Removal from Water Buffer, Neutral Sugars, and Certain Biological Solutions	231
3.0 ANALYSIS	
3.1 Western Blotting	235
3.1.1 Introduction	235
3.1.2 BioTrace™ NT Nitrocellulose Transfer Membrane	240
3.1.3 FluoroTrans® and BioTrace PVDF Membranes	248

TABLE OF CONTENTS

3.2 Affinity Activated or Activatable Membranes	259
3.2.1 Introduction	259
3.2.2 UltraBind™ Affinity Membrane	264
3.2.3 ELISA Assays Using Immunodyne® ABC Membranes	268
3.2.4 ELISA and Immunoassay Using AcroWell™ 96 Multi-Well Filter Plates with BioTrace PVDF and NT Membranes	272
3.2.5 ELISA and Immunoassay Using AcroWell 96 Multi-Well Filter Plates with Chemiluminescent (DELFI) Detection	282
3.2.6 Activation Protocol for Biodyne® C Membrane for Subsequent Covalent Linking of Ligands	292
4.0 PURIFICATION APPLICATIONS	
4.1 Mustang® Ion Exchange Membranes	297
4.1.1 Introduction	297
4.1.2 Purification on an Acrodisc® Unit with Mustang Q Membrane	298
4.1.3 Purification on an Acrodisc Unit with Mustang S Membrane	303
4.2 BioSeptra® Chromatography Resins	308
4.2.1 Introduction	308
4.2.2 Ceramic HyperD® Ion Exchange Resin	310
4.2.3 Blue Trisacryl® M Resin	325
4.2.4 IMAC HyperCel™ Resin for Immobilized Metal Ion Affinity Chromatography	332
4.2.5 HA Ultrogel® Resin	341
4.2.6 MEP HyperCel Resin for Purification of Antibodies by Hydrophobic Charge Induction Chromatography (HCIC)	350
4.2.7 Heparin HyperD Resin	359
4.2.8 Lysine HyperD Resin	365
4.2.9 Ultrogel AcA Size Exclusion Chromatography Resin	371
4.3 Rapid Purification Development/Scouting (Process Proteomics)	381
4.3.1 Introduction	381
4.3.2 Screening of Chromatographic Resin in an AcroPrep™ Multi-Well Plate Format	382
4.3.3 Column Scale-Up to Pilot Process Applications	390
4.4 Buffer Exchange, Desalting, and Concentration	391
4.4.1 Introduction	391
4.4.2 Microsep™ UF Spin Filter for Samples (1-3.5 mL)	393
4.4.3 Macrosep® UF Spin Filter for Samples (3.5-15 mL)	403
4.4.4 Jumbosep™ UF Spin Filter for Samples (15-60 mL)	413
4.5 Final Product Clarification	421
4.5.1 Introduction	421
4.5.2 Clarification of Samples (< 1 mL) by MF Filtration in Nanosep® Spin Filters	423
4.5.3 Filter Sterilization of Samples (1-100 mL) in Sterile Acrodisc Syringe Filters	426
5.0 SUPPORTING TECHNOLOGIES	
5.1 Water Supply	433
5.1.1 Introduction	433
5.1.2 Cascada™ BIO-Water System	434
5.1.3 Cascada AN-Water System	437
5.1.4 Cascada IX-Water System	440
5.1.5 Cascada LS-Water System	442
5.1.6 Cascada RO-Water System	444
5.2 Microbiology Testing	447
5.2.1 Monitoring for Bacteriological Contamination with MicroFunnel™ Plus Filter Funnels	447

6.0 TECHNICAL APPENDICES	
6.1 Evaluation of Chromatography Column Packing Efficiency	455
6.2 Optional Pre-Treatment to Improve Recovery of Samples from Ultrafiltration Spin Filters	458
6.3 Vacuum Manifold for Use with Multi-Well Filter Plates	459
6.4 General Filtration Procedures for AcroPrep™ and AcroWell™ Filter Plates	464
6.5 Chromatography Products Selection Guide	469
6.6 BioSeptra® Media	476
6.7 Protein Sample Preparation and Analysis Media	477
6.8 Chemical Compatibility Guide	480

TABLE OF CONTENTS

Notes: