

Subcellular Protein Fractionation Kit

78840

2131.0

Number	Description
78840	<p>Subcellular Protein Fractionation Kit, sufficient reagents for extracting 50 cell pellet fractions having packed cell volumes of 20 μl each (a total of ~2 g cell paste)</p> <p>Kit Contents:</p> <p>Cytoplasmic Extraction Buffer (CEB), 10 ml, store at -20°C</p> <p>Membrane Extraction Buffer (MEB), 10 ml, store at 4°C</p> <p>Nuclear Extraction Buffer (NEB), 10 ml, store at 4°C</p> <p>Pellet Extraction Buffer (PEB), 5 ml, store at room temperature</p> <p>Micrococcal Nuclease, ≥ 100 units/μl, 150 μl, store at -20°C</p> <p>Calcium Chloride (CaCl₂), 100 mM, 250 μl, store at 4°C</p> <p>Halt™ Protease Inhibitor Cocktail, 100X, 350 μl, store at 4°C</p> <p>Storage: Upon receipt store kit at -20°C, or store individual components as indicated above. Kit is shipped with dry ice.</p>

Introduction

The Thermo Scientific Subcellular Protein Fractionation Kit enables stepwise separation and preparation of cytoplasmic, membrane, nuclear soluble, chromatin-bound and cytoskeletal protein extracts from mammalian cultured cells or tissue. The first reagent added to a cell pellet causes selective cell membrane permeabilization, releasing soluble cytoplasmic contents. The second reagent dissolves plasma, mitochondria and ER/golgi membranes but does not solubilize nuclear membranes. After recovering the intact nuclei by centrifugation, a third reagent yields the soluble nuclear extract. A second nuclear extraction with micrococcal nuclease is performed to release chromatin-bound nuclear proteins. The recovered insoluble pellet is then extracted with the final reagent to isolate cytoskeletal proteins.

Extracts obtained with the Subcellular Protein Fractionation Kit are compatible with a variety of downstream applications including Western blotting, Thermo Scientific Pierce BCA Protein Assay (Product No. 23225), Thermo Scientific LightShift Chemiluminescent EMSA Kit (Product No. 20148), and reporter-gene and enzyme-activity assays. Extracts from each subcellular compartment generally have less than 15% contamination between fractions, which is sufficient purity for most experiments studying protein localization and redistribution.

Important Procedural Notes

- Thaw all buffers using a room temperature water bath, and keep CEB, MEB and NEB on ice until use. If precipitate occurs in PEB, mix vigorously to resuspend. Presence of a precipitate does not adversely affect PEB performance.
- Protease inhibitors are required to maintain extract integrity and function. Immediately before use, add protease inhibitors to CEB, MEB, NEB and PEB by diluting Halt Protease Inhibitor Cocktail 1:100 into each volume of buffer required.
- Perform all incubations at 4°C unless otherwise noted. Use a rotary shaker to avoid clumping of insoluble material during incubations.
- Perform all centrifugation steps at 4°C. Keep cell samples and extracts on ice unless otherwise noted.
- Subcellular protein extracts can be used directly in many downstream assays. Some applications might require dialysis or desalting to remove detergent and salts. Although the detergent in the MEB is not dialyzable, it does not interfere with isoelectric focusing. PEB contains a strong denaturing detergent that is not compatible with isoelectric focusing. For 2D analysis of cytoskeletal proteins, resuspend pellet directly in 2D sample buffer.

Additional Materials Required

- Ice-cold phosphate-buffered saline (PBS): 0.1 M sodium phosphate, 0.15 M sodium chloride; pH 7.2 (Product No. 28372)
- Rotary shaker in cold room
- Microcentrifuge at 4°C

Cell Culture Preparation

1. For adherent cells, harvest with trypsin-EDTA and then centrifuge at $500 \times g$ for 5 minutes. For suspension cells, harvest by centrifuging at $500 \times g$ for 5 minutes.
2. Wash cells by suspending the cell pellet with ice-cold PBS.
3. Transfer $1-10 \times 10^6$ cells to a 1.5 ml microcentrifuge tube and pellet by centrifugation at $500 \times g$ for 2-3 minutes.
4. Use a pipette to carefully remove and discard the supernatant, leaving the cell pellet as dry as possible.
5. Add ice-cold CEB containing protease inhibitors to the cell pellet (Table 1). Proceed to Subcellular Protein Extraction, using the reagent volumes indicated in Table 1.

Table 1. Reagent volumes for different packed cell volumes.*

<u>Packed Cell Volume (μl)</u>	<u>CEB (μl)</u>	<u>MEB (μl)</u>	<u>NEB (μl)</u>	<u>NEB (μl)</u> +CaCl ₂ , MNnase [†]	<u>PEB (μl)</u>
10	100	100	50	50	50
20	200	200	100	100	100
50	500	500	250	250	250
100	1,000	1,000	500	500	500

*For HeLa cells, 2×10^6 cells is equivalent to 20 μl packed cell volume. [†]MNase = Micrococcal Nuclease

Subcellular Protein Fractionation

Note: Scale this protocol depending on the cell pellet volume (Table 1). Maintain the volume ratio of CEB:MEB:NEB:PEB reagents at 200:200:100:100 μl, respectively.

1. After adding CEB to the cell pellet, incubate the tube at 4°C for 10 minutes with gentle mixing.
2. Centrifuge at $500 \times g$ for 5 minutes. Immediately transfer the supernatant (cytoplasmic extract) to a clean pre-chilled tube on ice.
3. Add ice-cold MEB containing protease inhibitors to the pellet. Vortex the tube for 5 seconds on the highest setting. Incubate tube at 4°C for 10 minutes with gentle mixing.
4. Centrifuge at $3,000 \times g$ for 5 minutes.
5. Transfer the supernatant (membrane extract) to a clean pre-chilled tube on ice.
6. Add ice-cold NEB containing protease inhibitors to the pellet. Vortex on the highest setting for 15 seconds. Incubate tube at 4°C for 30 minutes with gentle mixing.
7. Centrifuge at $5,000 \times g$ for 5 minutes. Transfer the supernatant (soluble nuclear extract) fraction to a clean pre-chilled tube on ice.
8. Prepare chromatin-bound extraction buffer by adding 5 μl of 100 mM CaCl₂ and 3 μl of Micrococcal Nuclease (300 units) per 100 μl of room temperature NEB.
9. Add room temperature NEB containing protease inhibitors, CaCl₂ and Micrococcal Nuclease to the pellet. Vortex on the highest setting for 15 seconds.
10. Incubate at room temperature for 15 minutes or in a 37°C water bath for 5 minutes.
11. After incubation, vortex on the highest setting for 15 seconds and centrifuge the tube at $16,000 \times g$ (highest setting of microcentrifuge) for 5 minutes.

12. Transfer the supernatant (chromatin-bound nuclear extract) fraction to a clean pre-chilled tube on ice.
13. Add room temperature PEB containing protease inhibitors to the pellet. Vortex on the highest setting for 15 seconds. Incubate at room temperature for 10 minutes.
14. Centrifuge the tube at $16,000 \times g$ (i.e., the highest microcentrifuge setting) for 5 minutes. Transfer the supernatant (i.e., the cytoskeletal extract) to a new tube.

Note: For same-day use, maintain fractions on ice for downstream applications and analysis. For long-term storage, store fractions at -80°C .

Troubleshooting

Problem	Possible Cause	Solution
Low cytoplasmic protein yield	Cells not lysed	Increase incubation time in CEB
	Cell pellet not dispersed	Mix cells gently during incubation
	CEB stored improperly	Store CEB at -20°C
Low membrane protein yield	Membranes solubilized with CEB	Decrease incubation time in CEB
	Incomplete membrane protein isolation	Increase time in MEB
Low soluble nuclear protein yield	Nuclei not extracted	Vortex thoroughly
	Incomplete nuclei isolation	Increase time of centrifugation after adding MEB
Low chromatin-bound protein yield	Calcium chloride or micrococcal nuclease not added	Add CaCl_2 and micrococcal nuclease to NEB before extraction
	Micrococcal nuclease stored improperly	Store micrococcal nuclease at -20°C
	Chromatin not completely degraded	Vortex thoroughly, add more micrococcal nuclease or incubate longer at 37°C
Low overall protein yield	Volumes of extraction reagents were not appropriate for given packed cell volume or tissue weight	Use the reagent volumes listed in Table 1
Proteins not compartmentalized	Incomplete lysis	Remove all PBS before adding CEB
		Vortex longer to completely disperse the pellet
		Increase incubation time
	Incomplete removal of extracts	Carefully remove all extract before proceeding to the next step
		Re-centrifuge sample and remove excess extract
		Rinse pellets with additional extraction buffers or PBS

Related Thermo Scientific Products

28372	BupH™ Phosphate Buffered Saline Pack, 40 packs
78439	Halt Protease Inhibitor Cocktail, EDTA-Free (100X), 10 ml
88216	Micrococcal Nuclease, ≥ 100 units/μl, 150 μl
78833	NE-PER Nuclear and Cytoplasmic Extraction Reagents, 1 kit
20148	LightShift® Chemiluminescent EMSA Kit
23224	Pierce BCA Protein Assay Reagent Kit
22660	Pierce 660 nm Protein Assay Reagent
89865	2-D Sample Preparation for Soluble Proteins
69558	Slide-A-Lyzer MINI Dialysis Units plus Float, 3.5K MWCO, 10 units
89882	Zeba™ Spin Desalting Columns, 0.5 ml, 25/pack

This product ("Product") is warranted to operate or perform substantially in conformance with published Product specifications in effect at the time of sale, as set forth in the Product documentation, specifications and/or accompanying package inserts ("Documentation") and to be free from defects in material and workmanship. Unless otherwise expressly authorized in writing, Products are supplied for research use only. No claim of suitability for use in applications regulated by FDA is made. The warranty provided herein is valid only when used by properly trained individuals. Unless otherwise stated in the Documentation, this warranty is limited to one year from date of shipment when the Product is subjected to normal, proper and intended usage. This warranty does not extend to anyone other than the original purchaser of the Product ("Buyer").

No other warranties, express or implied, are granted, including without limitation, implied warranties of merchantability, fitness for any particular purpose, or non infringement. Buyer's exclusive remedy for non-conforming Products during the warranty period is limited to replacement of or refund for the non-conforming Product(s).

There is no obligation to replace Products as the result of (i) accident, disaster or event of force majeure, (ii) misuse, fault or negligence of or by Buyer, (iii) use of the Products in a manner for which they were not designed, or (iv) improper storage and handling of the Products.

Current versions of product instructions are available at www.thermo.com/pierce. For a faxed copy, call 800-874-3723 or contact your local distributor.

© 2009 Thermo Fisher Scientific Inc. All rights reserved. Unless otherwise indicated, all trademarks are property of Thermo Fisher Scientific Inc. and its subsidiaries. Printed in the USA.