

Protocol: Identify Conditions that Solubilize Reversibly Aggregated Protein Samples


Material: aggregated protein sample (50 μ L for each protein sample to be tested) – up to six different protein samples can be applied with one kit.

Protocol for two aggregated protein samples:

1. Distribute 50 μ L aliquots from wells from OptiResc Reagent plate wells into large reservoir wells of OptiResc Inspection Plate (all 96 wells).
2. Transfer 2 x 0.5 μ L from each reservoir well of OptiResc Inspection Plate into small wells of OptiResc Inspection plate.
3. Add 0.5 μ L of aggregated protein sample A to all small wells (upper well a) and add sample B to all small wells (lower well b) within OptiResc Inspection Plate.
4. Seal OptiResc Inspection Plate with one Seal Sheet.
5. Inspect small wells with a microscope, i.e. a dissecting stereo microscope and note the wells that contain precipitate (aggregated protein solution) or clear drops (dissolved protein).
6. Employ OptiResc Reagent Listing (see reverse side of this brochure or www.dilyx.com/optiresc) to identify solutions that yield clear drops.

Scale up disaggregation reaction by mixing equal volumes of aggregated protein sample with solution that was identified with the OptiResc kit.

Consult the OptiResc Manual for details.

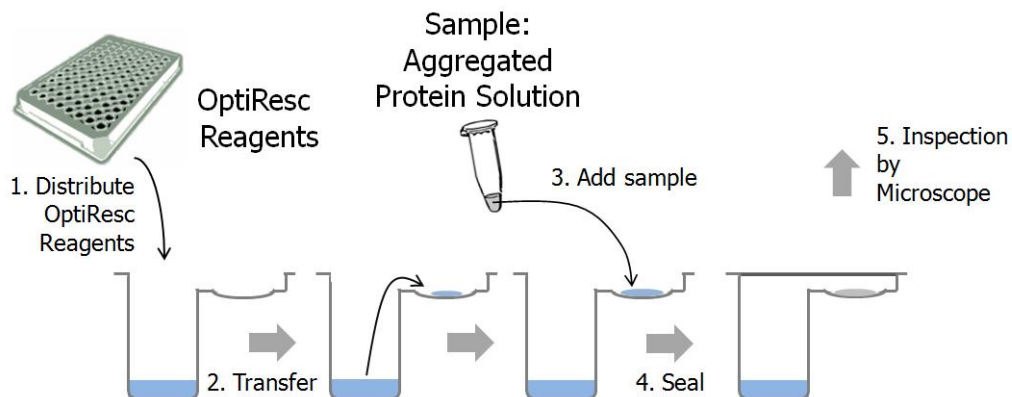

Product Information

Content:

- 1 x 96 well OptiResc Reagent Plate
- 1 x 96 x 2 well OptiResc Inspection Plate
- 3 x 96 well Seal sheets
- Instructions & Tech Sheet

Purpose

OptiResc Protein Aggregate Solubilization Kit

Systematic solution design and array-based deaggregation technology for:

- **Solubilizing reversibly aggregated Protein Samples**

For updated instructions and additional information please check: www.dilyx.com/optiresc

One OptiResc kit contains all consumable materials to identify solution conditions of up to six (6) different, reversibly aggregated protein samples.

Order Information

Order #: DLX-103-001

OptiResc Protein Aggregate Solubilization Kit

Price: \$ 599 (inquire about volume discounts)

Dilyx Biotechnologies, LLC
815 First Ave
Seattle WA 98104 USA

www.dilyx.com
info@dilyx.com
Phone: 425-205-0588
Fax: 202-207-0398

OptiResc Reagent Listing

Well		Buffer [#]			Additive		Well		Buffer [#]			Additive			
#	Row Col		Conc unit	pH	NAME	Conc unit	#	Row Col		Conc unit	pH	NAME	Conc unit		
1	A 1		Glycine	100 mM	3.0		49	E 1		Glycine	50 mM	3.0	Na ₂ SO ₄	500 mM	
2	A 2		Citric Acid	100 mM	3.2		50	E 2		Sodium Acetate	50 mM	4.5	Na ₂ SO ₄	500 mM	
3	A 3		PIPPS	100 mM	3.7		51	E 3		Bis-Tris	50 mM	6.0	Na ₂ SO ₄	500 mM	
4	A 4		Citric Acid	100 mM	4.0		52	E 4		MOPS	50 mM	7.0	Na ₂ SO ₄	500 mM	
5	A 5		Sodium Acetate	100 mM	4.5		53	E 5		Imidazole	50 mM	8.0	Na ₂ SO ₄	500 mM	
6	A 6		Na/K Phosphate	100 mM	5.0		54	E 6		CHES	50 mM	9.5	Na ₂ SO ₄	500 mM	
7	A 7		Sodium Citrate	100 mM	5.5		55	E 7		Citric Acid	50 mM	3.2	Arg/Glu*	50 mM	
8	A 8		Na/K Phosphate	100 mM	6.0		56	E 8		Na/K Phosphate	50 mM	5.0	Arg/Glu*	50 mM	
9	A 9		Bis-Tris	100 mM	6.0		57	E 9		ADA	50 mM	6.5	Arg/Glu*	50 mM	
10	A 10		MES	100 mM	6.2		58	E 10		HEPES	50 mM	7.5	Arg/Glu*	50 mM	
11	A 11		ADA	100 mM	6.5		59	E 11		Tris	50 mM	8.5	Arg/Glu*	50 mM	
12	A 12		Bis-Tris Propane	100 mM	6.5		60	E 12		CAPS	50 mM	10.0	Arg/Glu*	50 mM	
13	B 1		Ammonium Acetate	100 mM	7.0		61	F 1		Glycine	50 mM	3.0	Tween 20	1 % (w/v)	
14	B 2		MOPS	100 mM	7.0		62	F 2		Sodium Acetate	50 mM	4.5	Tween 20	1 % (w/v)	
15	B 3		Na/K Phosphate	100 mM	7.0		63	F 3		Bis-Tris	50 mM	6.0	Tween 20	1 % (w/v)	
16	B 4		HEPES	100 mM	7.5		64	F 4		MOPS	50 mM	7.0	Tween 20	1 % (w/v)	
17	B 5		Tris	100 mM	7.5		65	F 5		Imidazole	50 mM	8.0	Tween 20	1 % (w/v)	
18	B 6		EPPS	100 mM	8.0		66	F 6		CHES	50 mM	9.5	Tween 20	1 % (w/v)	
19	B 7		Imidazole	100 mM	8.0		67	F 7		Citric Acid	50 mM	3.2	Solubilisin TM	100 % (w/v)	
20	B 8		Bicine	100 mM	8.5		68	F 8		Na/K Phosphate	50 mM	5.0	Solubilisin TM	100 % (w/v)	
21	B 9		Tris	100 mM	8.5		69	F 9		ADA	50 mM	6.5	Solubilisin TM	100 % (w/v)	
22	B 10		CHES	100 mM	9.0		70	F 10		HEPES	50 mM	7.5	Solubilisin TM	100 % (w/v)	
23	B 11		CHES	100 mM	9.5		71	F 11		Tris	50 mM	8.5	Solubilisin TM	100 % (w/v)	
24	B 12		CAPS	100 mM	10.0		72	F 12		CAPS	50 mM	10.0	Solubilisin TM	100 % (w/v)	
25	C 1		Glycine	50 mM	3.0	NaCl	150 mM	73	G 1		Glycine	50 mM	3.0	Glycerol	20 % (w/v)
26	C 2		Sodium Acetate	50 mM	4.5	NaCl	150 mM	74	G 2		Sodium Acetate	50 mM	4.5	Glycerol	20 % (w/v)
27	C 3		Bis-Tris	50 mM	6.0	NaCl	150 mM	75	G 3		Bis-Tris	50 mM	6.0	Glycerol	20 % (w/v)
28	C 4		MOPS	50 mM	7.0	NaCl	150 mM	76	G 4		MOPS	50 mM	7.0	Glycerol	20 % (w/v)
29	C 5		Imidazole	50 mM	8.0	NaCl	150 mM	77	G 5		Imidazole	50 mM	8.0	Glycerol	20 % (w/v)
30	C 6		CHES	50 mM	9.5	NaCl	150 mM	78	G 6		CHES	50 mM	9.5	Glycerol	20 % (w/v)
31	C 7		Citric Acid	50 mM	3.2	NaCl	500 mM	79	G 7		Citric Acid	50 mM	3.2	Betaine	2 M
32	C 8		Na/K Phosphate	50 mM	5.0	NaCl	500 mM	80	G 8		Na/K Phosphate	50 mM	5.0	Betaine	2 M
33	C 9		ADA	50 mM	6.5	NaCl	500 mM	81	G 9		ADA	50 mM	6.5	Betaine	2 M
34	C 10		HEPES	50 mM	7.5	NaCl	500 mM	82	G 10		HEPES	50 mM	7.5	Betaine	2 M
35	C 11		Tris	50 mM	8.5	NaCl	500 mM	83	G 11		Tris	50 mM	8.5	Betaine	2 M
36	C 12		CAPS	50 mM	10.0	NaCl	500 mM	84	G 12		CAPS	50 mM	10.0	Betaine	2 M
37	D 1		Glycine	50 mM	3.0	Trehalose	1.0 M	85	H 1		H2O	100 %			
38	D 2		Sodium Acetate	50 mM	4.5	Trehalose	1.0 M	86	H 2		H2O	100 %			
39	D 3		Bis-Tris	50 mM	6.0	Trehalose	1.0 M	87	H 3						
40	D 4		MOPS	50 mM	7.0	Trehalose	1.0 M	88	H 4					AmSulfate	3 M
41	D 5		Imidazole	50 mM	8.0	Trehalose	1.0 M	89	H 5					Acetonitrile	80 % (v/v)
42	D 6		CHES	50 mM	9.5	Trehalose	1.0 M	90	H 6		PEG 1450	10 %		NaCl	50 mM
43	D 7		Citric Acid	50 mM	3.2	TMAO	500 mM	91	H 7					DDT	1 mM
44	D 8		Na/K Phosphate	50 mM	5.0	TMAO	500 mM	92	H 8					DDT	5 mM
45	D 9		ADA	50 mM	6.5	TMAO	500 mM	93	H 9					DDT	15 mM
46	D 10		HEPES	50 mM	7.5	TMAO	500 mM	94	H 10					BME	2.5 mM
47	D 11		Tris	50 mM	8.5	TMAO	500 mM	95	H 11					BME	10 mM
48	D 12		CAPS	50 mM	10.0	TMAO	500 mM	96	H 12					BME	20 mM

TMAO, Trimethylamine N-Oxide; PIPPS, Piperazine-N, n'-Bis (3-Propanesulfonic Acid); MES, 2-(N-morpholino) ethanesulfonic acid; MOPS, 3-(N-morpholino) propanesulfonic acid; HEPES, 4-(2-hydroxyethyl)-1-piperazineethanesulfonic acid; Arg/Glu*: 50mM of each Arginine and Glutamate; DDT, DL-Dithiothreitol; BME, 2-Mercaptoethanol; Betaine, Trimethyl-Glycine; CAPS, N-cyclohexyl-3-amino-propanesulfonic acid; ADA, N-(2-Acetamidimidino)diacetic Acid; Tris, tris(hydroxymethyl)aminomethane; CHES, 2-(N-Cyclohexylamino)ethane Sulfonic Acid; EPPS, N-(2-hydroxyethyl)piperazine-N'-(3-propanesulfonic acid). # pH values for buffers used only; * each amino acid is 50 mM © Dilyx Biotechnologies, LLC; Patent pending