

SuperSpin™ Desalator

Data and Instructions



Based on the principles behind size exclusion chromatography, cross-linked neutral polysaccharide particles with very small pores are packed into spin tubes for rapid desalting and / or buffer exchange. SuperSpin Desalator achieves desalting or buffer exchange in just a few minutes. Up to 24 samples (depending on the type of microcentrifuge) of 10 – 100 µl can be processed in one spin. It is a much faster and more efficient approach when compared to dialysis tubes or membrane ultrafiltration. The small porous particles provide huge

surface area with very short diffusion distance, which means small molecules such as salt can be partitioned rapidly. In comparison, both dialysis tube and membrane have very low surface area. It always takes much longer to conduct dialysis (typically from a few hours to a few days). Membrane ultrafiltration of small samples always experiences membrane blockage and severe loss of valuable materials.

Key benefits:



- Rapid desalting or buffer exchange
- Very little loss of target molecules (typically > 95% recovery)
- Most proteins (> 6,000 dalton) and DNAs (> 10 bp) can be desalted
- DNase free

SuperSpin desalator is particularly useful for the following applications:

- Desalting of histidine-tagged proteins (e.g. imidazole and NaCl) recovered from IMAC SuperSpin
- Desalting of samples before loading to SDS-PAGE
- Desalting of samples before conducting other analysis
- Buffer exchange, for example, after low pH elution
- Desalting of DNAs

Operational instruction

1. Place a SuperSpin Desalator into a standard 1.5 ml microcentrifuge tube. (Caution: don't open the lid of SuperSpin Desalator in this stage). Keep the cap of the 1.5 ml tube open, or cut it off if it interferes the spin process.
2. Spin at 6500 rpm for 10 seconds to pack down the particles.
3. Gently open the lid of the spin tube. Slowly load 200 µl of the equilibration buffer of choice. Close the lid and spin at 6500 rpm for 30 seconds. *The liquid retention time in the spin tube may vary for different buffers. The resin bed needs be fully dried in this spin step. By visual checking, the resin bed should become white dry. Otherwise, spin for another 30 seconds.*
4. Empty the 1.5 ml centrifuge tube and replace in the Desalator tube. Repeat the above Step 3 one more time.
5. Place the spin tube into a fresh 1.5 ml tube. Load the sample upon top of the particles gently. The loading volume is 10 µl to 100 µl. Close the spin tube lid and spin at 6500 rpm for 30 seconds.
6. The liquid collected in the 1.5 ml tube is the final desalted sample.

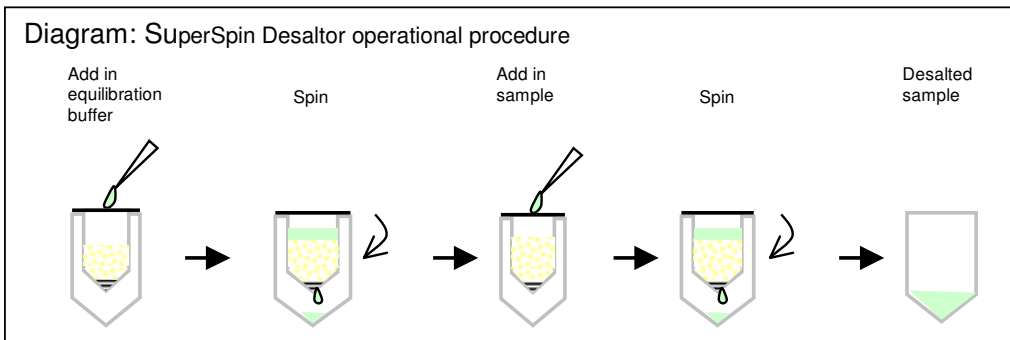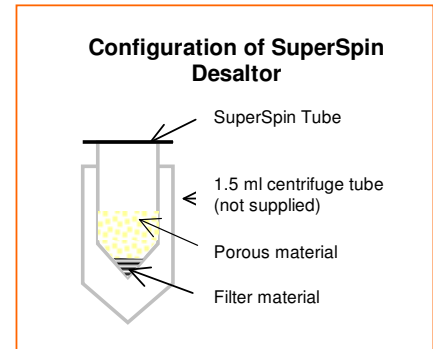

BIOTOOLOMICS

Technical data

Biological samples with molecule size > 6,000 dalton can be desalted with SuperSpin Desaltor.

Protein	Sample loading	Salt removal	Protein recovery
Lysozyme (14.6K, 1 mg/ml in 10 mM Tris/HCl plus 1 M NaCl)	10 µl	100%	99%
Lysozyme (14.6K, 1 mg/ml in 10 mM Tris/HCl plus 1 M NaCl)	100 µl	86.5%	96%
BSA (66K, 1 mg/ml in 10 mM Tris/HCl plus 1 M NaCl)	10 µl	100%	83%
BSA (66K, 1 mg/ml in 10 mM Tris/HCl plus 1 M NaCl)	100 µl	83.5%	98%

Ordering information

Product	Quantity	Code no.
SuperSpin Desaltor	50	210101

Related products	Quantity	Code no.
Ni SuperSpin	50	150101
Cu SuperSpin	50	150102
Co SuperSpin	50	150103
Zn SuperSpin	50	150104

BioToolomics Limited
Consett Innovation Centre
Ponds Court Business Park
Genesis Way
Consett
County Durham DH8 5XP
UK

www.biotooolomics.com

Registered or registration-pending trademark of BioToolomics Ltd: BioToolomics, SepFast and SuperSpin. Triton is a registered trademark of Union Carbide inc. Tween is a registered trademark of ICI Americas Inc.

All goods and services are sold subject to the terms and conditions of sale of BioToolomics Ltd. The company reserves the rights, subject to regulatory or contractual approval, if required, to make changes in the specifications and features shown herein, or discontinue the products described at any time without notice or obligation. Contact BioToolomics Ltd for the most current information.

© 2006-2008 BioToolomics Ltd – All rights reserved.